


Installation of EP3View 2.35 with EP4 Model

This is a non free software version. First time you run the software you need to contact Nanofilm to get a software key.

Minimum system requirements

- 2.0 GHz Pentium processor or equivalent
- 1GB RAM
- 32-bit true-color graphics adapter, 1280 × 1024 pixels
- CD-ROM drive
- 2 USB interfaces (dongle, joystick)
- 2 × 100 MBit Fast Ethernet network adapters (EP3, corporate LAN)
- Windows 2000® (Service Pack 3 or higher) or Windows XP® or Vista
- Microsoft DirectX 8.1 for Logitech "Wingman" joystick (*optional*)

- 1) Run the installer and follow the online instructions. Destination folder must be "C:\Accurion".
- 2) When you get prompted to restart the computer select "restart later"!!
- 3) Follow the online instructions to install all software components
- 4) Copy the latest user from "D:\Backup\EP3View V2" or another source to "C:\Accurion\EP3View V2"
- 5) Restart the system.
- 6) Find the name of your computer by right click on "my computer" on your desktop and select properties. Copy the name under the tab "computer name\full computer name".
- 7) Open the file "C:\Accurion\EP3View V2\tools\EP4Model via Viserver\EP3ToEP4.ini" with notepad and paste your computer name as host (see marking).




```

[general]
UseEP4Model=True
//if no dongle, disable EP3 Model usage (this VI only)
[EP4Model]
Host="pcep3"
Port=3364
ReceivingVI="mod_CrossApp.vi"

```

8) Open the file “C:\Accurion\EP4Model\program\model\model.ini.xml” with notepad and paste your computer name as host (see marking):

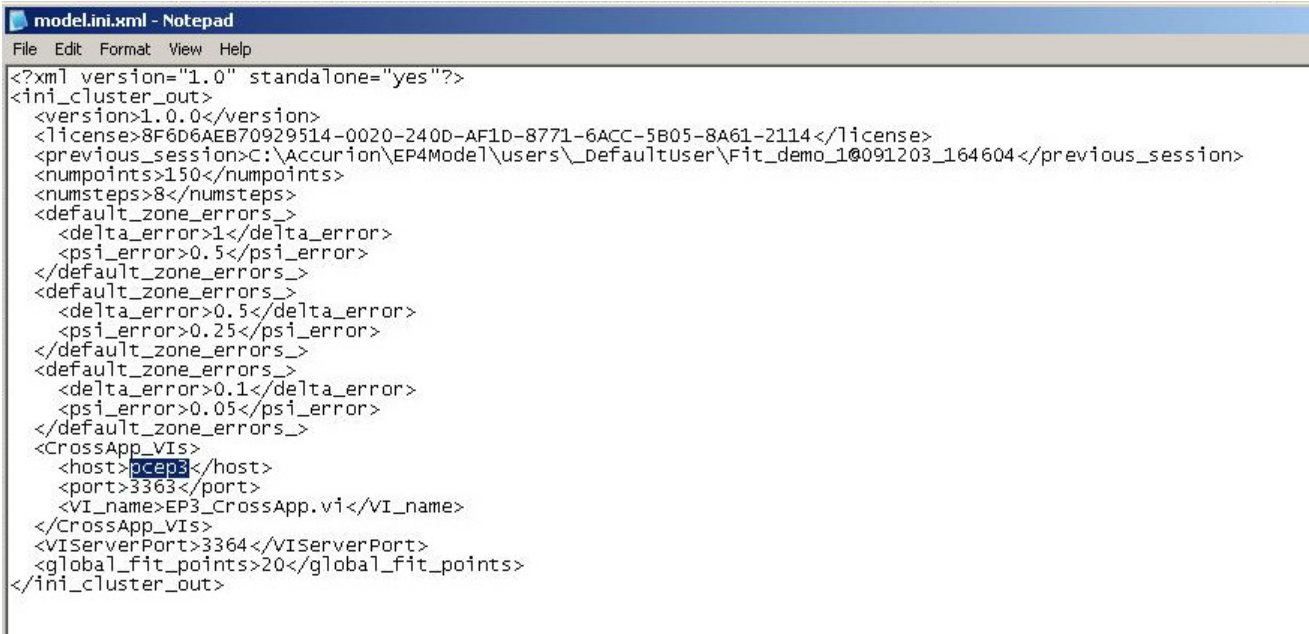

```

model.ini.xml - Notepad
File Edit Format View Help
<?xml version="1.0" standalone="yes"?>
<ini_cluster_out>
  <version>1.0.0</version>
  <license>8F6D6AEB70929514-0020-240D-AF1D-8771-6ACC-5B05-8A61-2114</license>
  <previous_session>C:\Accurion\EP4Model\users\_Defaultuser\Fit_demo_1@091203_164604</previous_session>
  <numpoints>150</numpoints>
  <numsteps>8</numsteps>
  <default_zone_errors_>
    <delta_error>1</delta_error>
    <psi_error>0.5</psi_error>
  </default_zone_errors_>
  <default_zone_errors_>
    <delta_error>0.5</delta_error>
    <psi_error>0.25</psi_error>
  </default_zone_errors_>
  <default_zone_errors_>
    <delta_error>0.1</delta_error>
    <psi_error>0.05</psi_error>
  </default_zone_errors_>
  <CrossApp_VIS>
    <host>accp3</host>
    <port>3363</port>
    <VI_name>EP3_CrossApp.vi</VI_name>
  </CrossApp_VIS>
  <VIsServerPort>3364</VIsServerPort>
  <global_fit_points>20</global_fit_points>
</ini_cluster_out>

```

9) Start EP3View and EP4Model, copy the license request code you get and contact Accurion to get a new licence key.